


Ivium MCF Cell

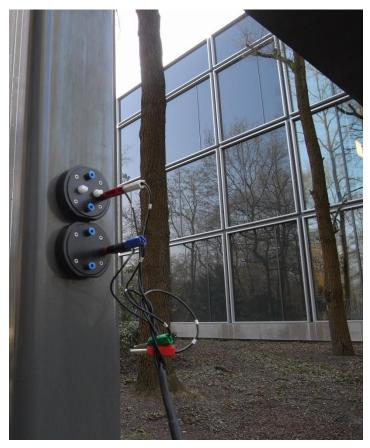

· In the Field In the Laboratory

Coating resistance: EIS measurements without damaging the coating!

The MCF cell is magnetic and can be clamped to objects in any position. In this way the object constitutes the bottom of the flat-cell, and the object itself can be used as electrode. The cell can be filled using a tube that leads directly into the cell. A similar tube allows air to escape to ensure a completely filled cell. A rubber seal prevents electrolyte leakage. The back of the cell consists of a stainless steel counter electrode. If desired a reference electrode can be placed near the electrode using either one of 2 holes designed for this purpose.

Multi-Functional Magnetic Flat cell:

- Magnetic
- Electrode
- Flat cell
- Can be clamped:

to any steel object in any position

- Applications:
 - Corrosion
 - Coatings
 - Materials






THE NETHERLANDS:

Ivium Technologies B.V. De Zaale 11 5612 AJ Eindhoven The Netherlands

tel. +31 40 2390600 fax. +31 40 2390601 e-mail info@ivium.eu www.ivium.com

Ivium Technologies USA 961687 Gateway Blvd., Suite 201D Fernandina Beach, FL 32034


phone: 800-303-3885 (toll free) / 904-310-9060 (office) fax: 904-310-9068 e-mail pete@ivium.us www.ivium.us

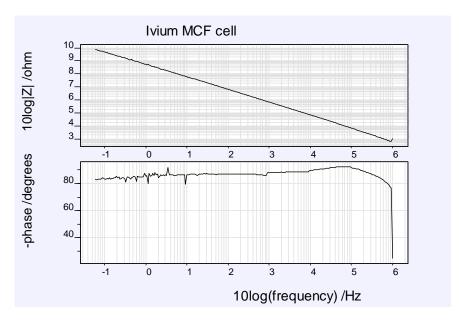

EIS measurement of coated Stainless Steel 416; CompactStat and Ivium MCF cell, 1MHz-0.06Hz, amplitude 20mV, rainwater electrolyte

Specifications

Bottom view Top view

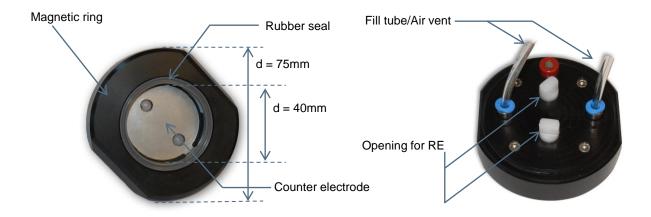

Dimensions

Size MCF cell

outer diameter: 75mm Fill tube: 4 mm/2.2mm height: 30mm RE opening: 6.2mm

Cell compartment

diameter: 40mm height: 10mm volume: 12.5 cm³

Weight 150 gram

Materials Cell – POM Electrode – Steel 316L

Seals – Silicone rubber